

# Technical Data Sheet:

## CULTEX® RFS

### Radial Flow System for Cell-based Exposure



*The Airborne Exposure Experts*



This is a high-end exposure system for the direct exposure of cell culture inserts or Petri dishes. The modular system consists of two main parts, the aerosol-guiding module and the sampling module housing the cell culture inserts, or Petri dishes.

The aerosol-guiding module can be fitted tightly to the sampling module, thus realizing a close connection between the two parts before exposure. The sampling module houses three cell culture inserts (or Petri dishes), which can be separately supplied with medium.

## Exposure systems

### General features

- Direct exposure technology at the air-liquid interface of cell cultures
- Cellular systems: cell lines and primary cells from the respiratory tract as mono- or co-cultures
- Test compounds
  - Gases
  - Particles (fine to nanoparticles)
  - Complex mixtures (e.g. tobacco smoke)
- Application areas
  - Industrial chemicals (particulate & gaseous compounds)
  - Consumer products
  - Tobacco smoke
  - E-cigarette vapor
- Pharmaceutical and therapeutic products
- Pesticides
- In- and outdoor analysis
- Clients
  - Universities
  - Regulatory Bodies
  - Military
  - Pharmaceutical, chemical and tobacco industry
  - Contract research laboratories

## Basic principles of the exposure systems

### Module design

- Inlet adapter
- Exposure top – Aerosol guiding module
- Base module – Sampling module
- Socket module
- Locking module

### Dilution

- Use of the system with or without dilution

### Test atmosphere

- One sampling point
- Radial flow system
- Homogenous distribution of the test atmosphere to the surface of the cells
- Reproducible deposition of the particles on the cell surface
- Establishment of dose-response relationships
- Repeated exposure studies

### Biological test system

#### Cells

- Maintenance of cell viability
  - 37°C by a temperature-controlled flow of water
  - Medium supply from below the insert membrane (static, intermittent, continuous)
- Application of 6.5 mm, 12 mm and 24 mm Falcon® or Transwell® inserts via special adapters

#### Bacteria


- Special adapter for 35 mm Petri dishes for running a modified Ames assay

# Technical Data Sheet: CULTEX® RFS – Radial Flow System for Cell-based Exposure

## Air-liquid interphase – *in vitro* exposure

### Exposure systems – CULTEX® RFS – Radial Flow System

- Flexible design of the exposure top with aerosol inlets for commercial available membrane inserts (Falcon®/Transwell®) of different size and 35 mm Petri dishes
- 3 exposure positions
- Individual exposure chambers (3 membrane inserts)
- Tempering of the exposure chambers
- Individual medium supply of the independent exposure chambers

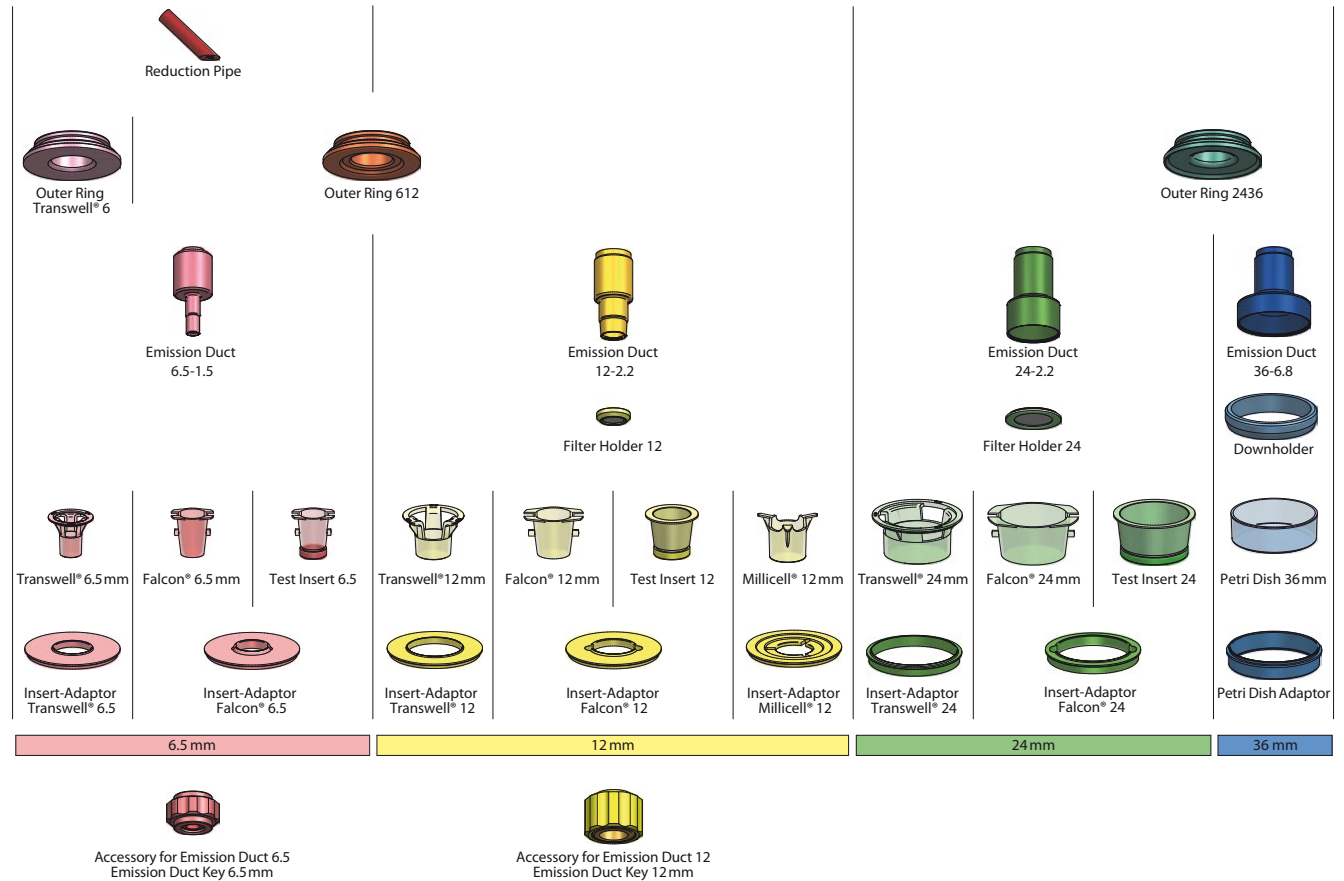

The Airborne Exposure Experts

**Contact us:**  
**Cultex® Laboratories GmbH** · Feodor-Lynen-Straße 21 · 30625 Hannover · Germany · Phone: +49 (0)511 563586-0 · [www.cultex-laboratories.com](http://www.cultex-laboratories.com) · [info@cultex-laboratories.com](mailto:info@cultex-laboratories.com)  
 Our products are subject to continuous development and we reserve the right to amend specifications without notice.  
 © 2016 Cultex® Laboratories GmbH. All rights reserved. Version 1.1 | Design by ahlersheinel.com